




Illumina Proactive connectivity guide

Easily connect your MiSeq™ i100 Series Sequencing System

Connecting your MiSeq i100 Series sequencing system ensures smoother operation, enhanced efficiency, and prompt resolution when support is needed.

Follow these steps to get connected

- From the **Start Menu** (1) on the Home Screen, select **Settings**, then select **Cloud settings** (2)
- Under the **Illumina Proactive** menu (3), select the checkmark to send instrument performance data to Illumina
- Click **Save**

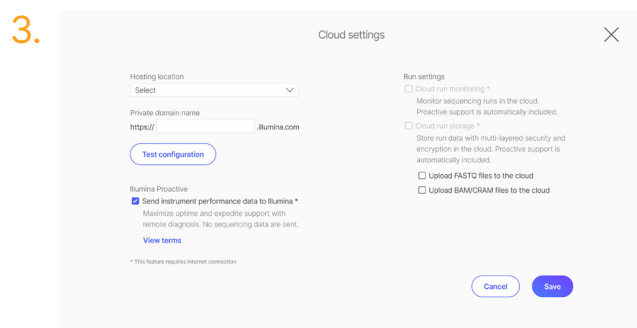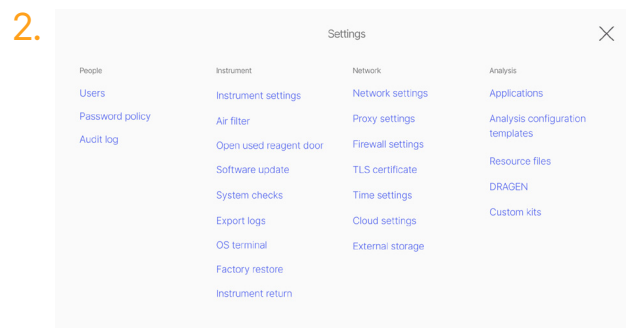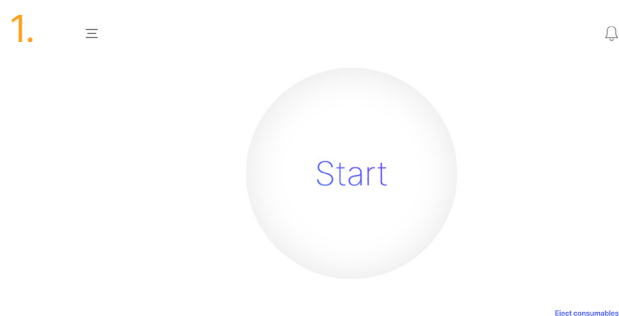

For more information, contact your local field applications scientist or field service engineer

Illumina • 1.800.809.4566 toll-free (US) • +1.858.202.4566 tel • techsupport@illumina.com • www.illumina.com

© 2025 Illumina, Inc. All rights reserved. All trademarks are the property of Illumina, Inc. or their respective owners. For specific trademark information, see www.illumina.com/company/legal.html. M-GL-03317 v1.0